## Instructions for using the Switch kit with the Timing ramp

The 3841 and 2341 Switch kits are designed for use with the Timing ramp (3842). A kit contains:

- Two magnetic proximity switches that fit into the holes in the ramp.
- A magnet which attaches to the car using the sticky pads supplied.

The **3841** Switch kit has mini din plugs for use with a data logger that will accept two digital Smart Q sensors e.g. EasySense Q3+, VISION, V-Log or V-Hub.


Note: As two inputs are required this kit is **not** suitable for use with EasySense 1Link.


The **2341** Switch kit has jack plugs and is suitable for use with the Vu or Vu+ logger.


The switches will detect the magnet /car as it travels down the ramp.

Assemble the Timing ramp according to its instruction leaflet.




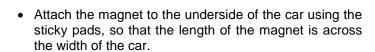

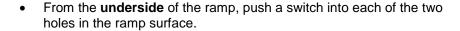

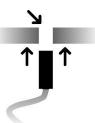


Make sure the switch is flush with the upper surface of the ramp

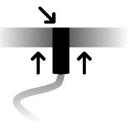




- Connect the switch that will be used first for timing to the input labelled A on the data logger.
- Connect the switch used second to input B on the data logger.

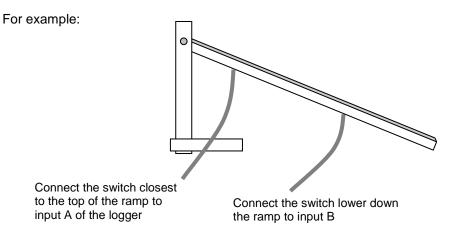

In the EasySense software (or on the data logger), select Timing and then how you want to record the data. You can use the switches to measure:

- The **Time** the car took to travel from switch A to switch B.
- The **Speed** of the car travelling between switch A and switch B (enter the distance between the two switches as the parameter e.g. 0.5 m).
- The car's **Acceleration** as it travels from A to B (enter the length of the magnet that will pass over the switches as the parameter e.g. 19 mm).

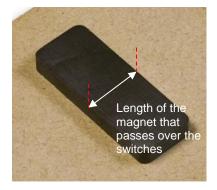

## **Limited warranty**

For information about the terms of the product warranty, see the Data Harvest website at: https://data-harvest.co.uk/warranty.



WEEE (Waste Electrical and Electronic Equipment) Legislation

Data Harvest Group Ltd is fully compliant with WEEE legislation and is pleased to provide a disposal service for any of our products when their life expires. Simply return them to us clearly identified as 'life expired' and we will dispose of them for you.